

## WB Developing and Fixing Kit

**Catalog#** BWR1070

**Size:** 100 ml (RTU)

**Lot #** Check on the product label

### Introduction

WB Developing and Fixing Kit is a processing kit, which can be used for Western, Northern and Southern Blotting after squashed. This kit is especially suitable for X-ray developing and fixing, as well as normal black & white negative film or photograph. This kit can be used solely without any other assistant reagent.

### Kit Components

Components	Size	Storage Instruction
Developing Solution	100 ml	Store at room temperature in dark for 3 months, or at 4°C in dark for long term.
Fixing Solution	100 ml	Store at room temperature for 3 months, or at 4°C for long term.

### Protocol

1. Developing: Immerse X-ray film in the Developing Solution for 10 seconds to 10 minutes. End user can grasp the time by observing under red light, stop developing once it achieves the experimental purpose.
2. Wash: Wash film with clean water to remove Developing Solution completely.
3. Fixing: Immerse the film in the Fixing Solution for 3-5 min. No data indicate that long fixing time affects the efficiency of film processing.
4. Wash: Wash film with clean water to remove various solutions thoroughly.

FOR RESEARCH USE ONLY, NOT FOR DIAGNOSTIC AND CLINICAL USE.

Chongqing Biospes Co., Ltd Tel: +86-23-67567091 Fax: +86-23-67745923

7F, Bldg B, High-tech Venture Park, # 107 Erlang Chuangye Rd, Jiulongpo District, Chongqing, 400039, China

[www.biospes.com](http://www.biospes.com)